


SMART SHUTTER INSTALLATION


STEP 01

REMOVE FILTER WHEEL COVER PLATE

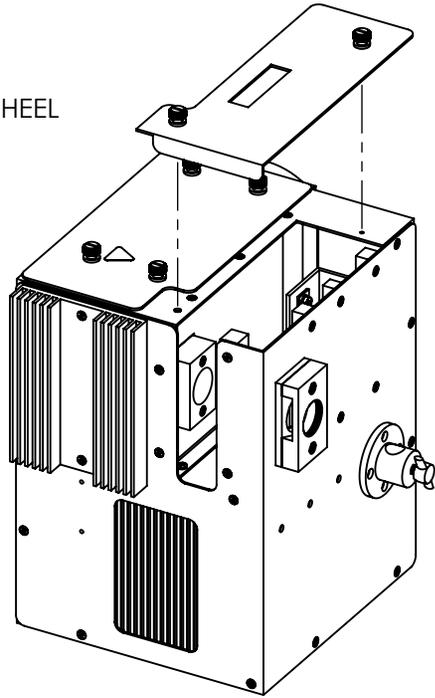


STEP 02

REMOVE DROP-IN FILTER HOLDER BY REMOVING THE 2 MOUNTING SCREWS

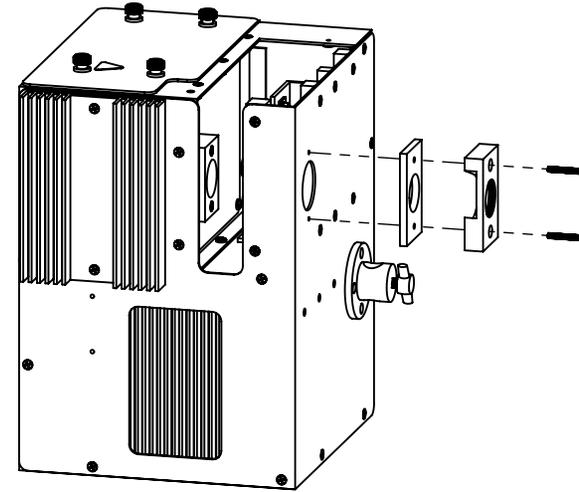


STEP 03

ATTACH MALE TO MALE C-MOUNT TO SMART-SHUTTER WITH THE DOVETAIL PORTION ON THE OUTSIDE.

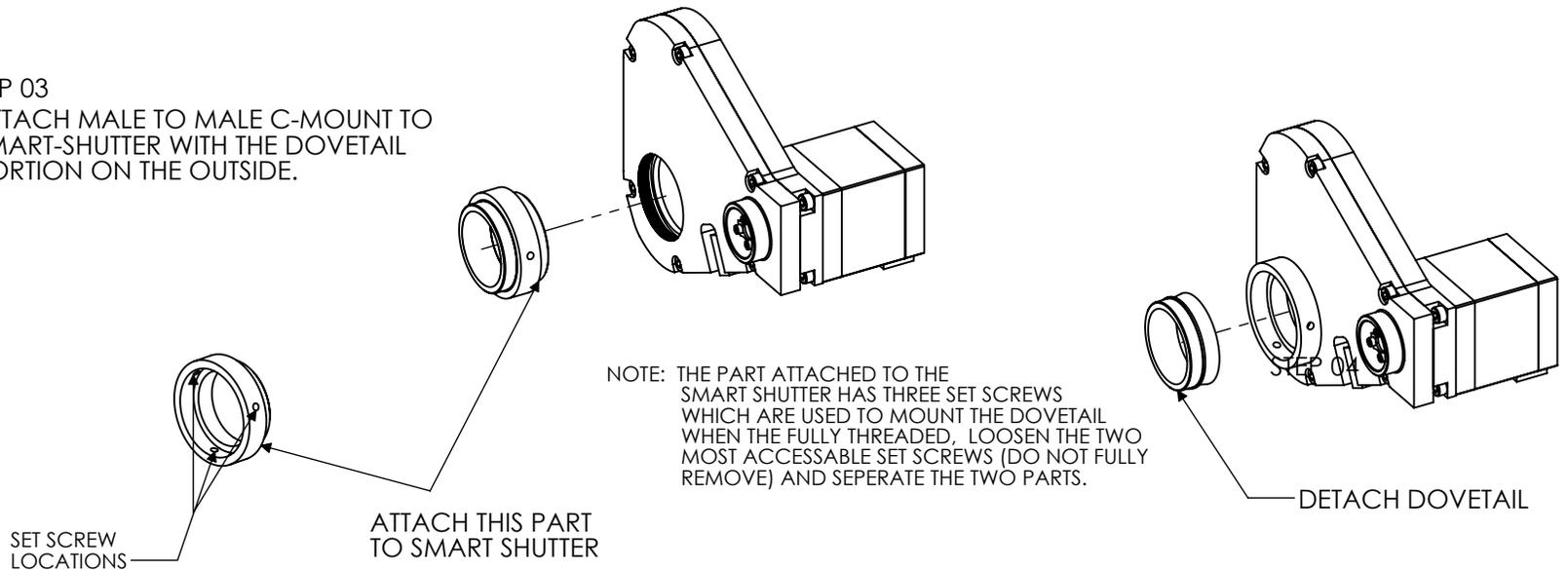


SMART SHUTTER INSTALLATION


STEP 05
ATTACH DOVETAIL TO
C-MOUNT BLOCK INSIDE
THE CABINET

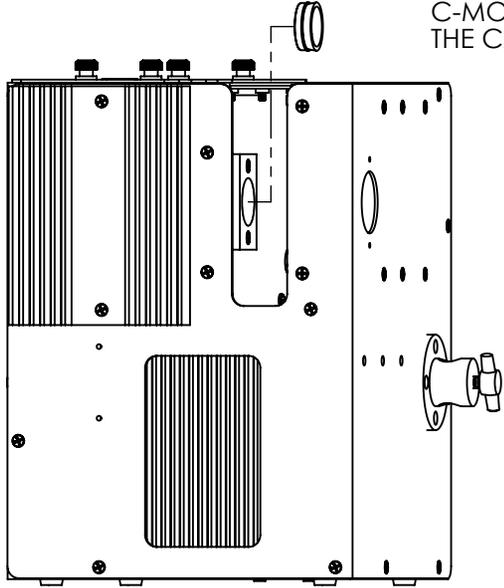


STEP 06
INSERT SMART SHUTTER SO THE TWO PARTS
OF THE C-MOUNT ADAPTER ARE ALIGNED.

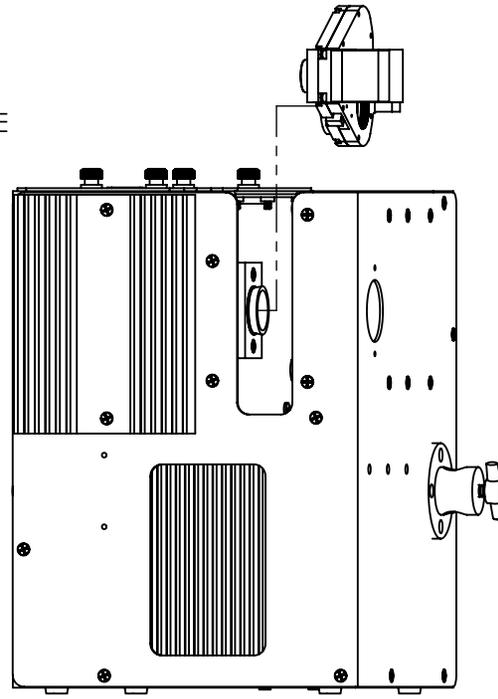

STEP 07
TIGHTEN THE TWO
EXPOSED SET SCREWS
TO MOUNT THE SMART SHUTTER

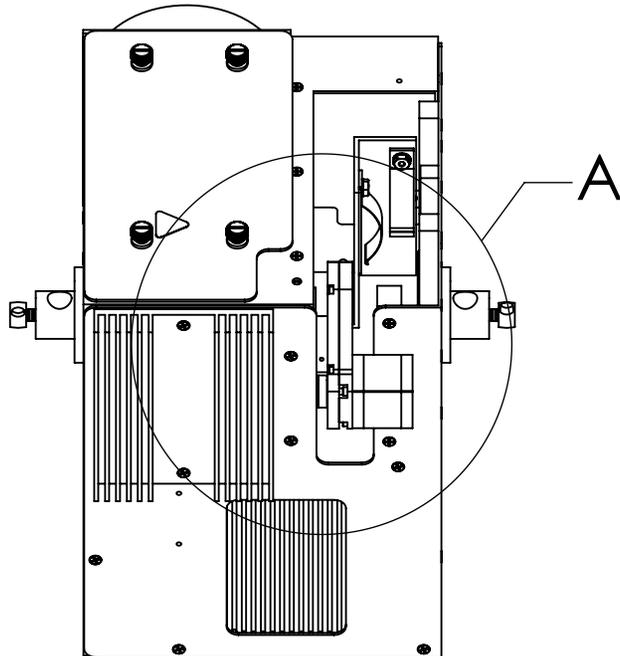

DETAIL A
SCALE 1 : 2

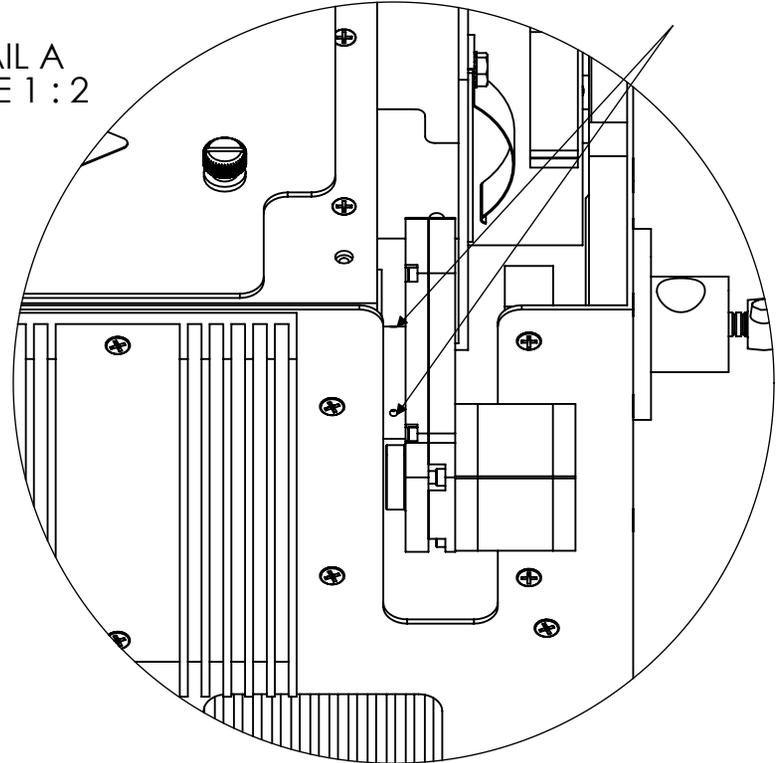

SMART SHUTTER INSTALLATION

